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Instructions for Use of TPP Vacuum Filtration «rapid»- Filtermax

The TPP filtration system «rapid»-Filtermax is suitable for sterile filtration of cell culture media, sera and aqueous solution.

The 0.22 µm polyether sulfone (PES) membrane ensures high flow rates with low protein binding and foam formation. The «rapid»-Filtermax is available as a filtration unit (filter top incl. reservoir flask) as well as a separate filter top or flask, with a DIN GL45 thread.

The filtration system is available in volume sizes 150, 250, 500 and 1000 ml. The individual system components are all identical in construction, available separately and can be combined with each other.

The square foot of the reservoir ensures a safe stand during sterile filtration.

The «rapid»-Filtermax is intended for single use only.

Note:

Follow national regulations when handling biological material, use appropriate protective clothing

Take note of the regulations of aseptic working during the work process.

Instructions:

- Check the packaging material and the product for intactness. Use only faultless material.
- Check the expiry date (EXP) on label and packaging. Only use products with valid EXP.
- Open the packaging in a sterile environment and remove a complete system for use. Only

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Advice and Recommendations:

- *The graduation serves as a reference point for the filling quantity and is not an absolute value.*
- *Alcohol 90 % with mechanical stress (rubbing) can dissolve or blur the white colour of the lettering.*

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Information

- The user should provide following materials:
Vacuum source; vacuum tubing.
- If a solution contains heavy particles to be filtered, it perform a pre-filtering using a fiberglass system.
- Solutions with a low particle density and a high filtration throughput can be attached to another reservoir for additional filtration. Please observe the rules of aseptic work.
- When filtering very dilute solutions, binding studies should be performed beforehand.
- Once the pouch is open, keep the product in a sterile area to ensure sterility.
- The flask is not suitable for freezing use.
- Do not autoclave product.
- Do not use the Filtermax in direct patient care applications or diagnostic procedures; it is intended for research use only

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To avoid possible injury due to implosions:

- Do not exceed 700 mm Hg differential vacuum at 25 °C
- Filter top: Only use sterile plastic or glass bottles that are approved for sterile filtration under vacuum. These should be free of any damage (danger of implosion). Pay attention to the max. total pressure.
- Wear eye-protection and perform the filtration in special safety units.
- Do not use hand-held vessels while filtering under vacuum.

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Technical Data:

Materials

Filter-membrane	Hydrophilic polyether sulfone (PES), type TPP fast flow
Flask and top	PS
Screw cap	PE
Vacuum connector	PP

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Additional:

Instructions for use, chemical resistance lists, and quality certificates of the respective products are available for download on the homepage www.tpp.ch